


Twist Bottle Adapter

Part Number PUA1507T

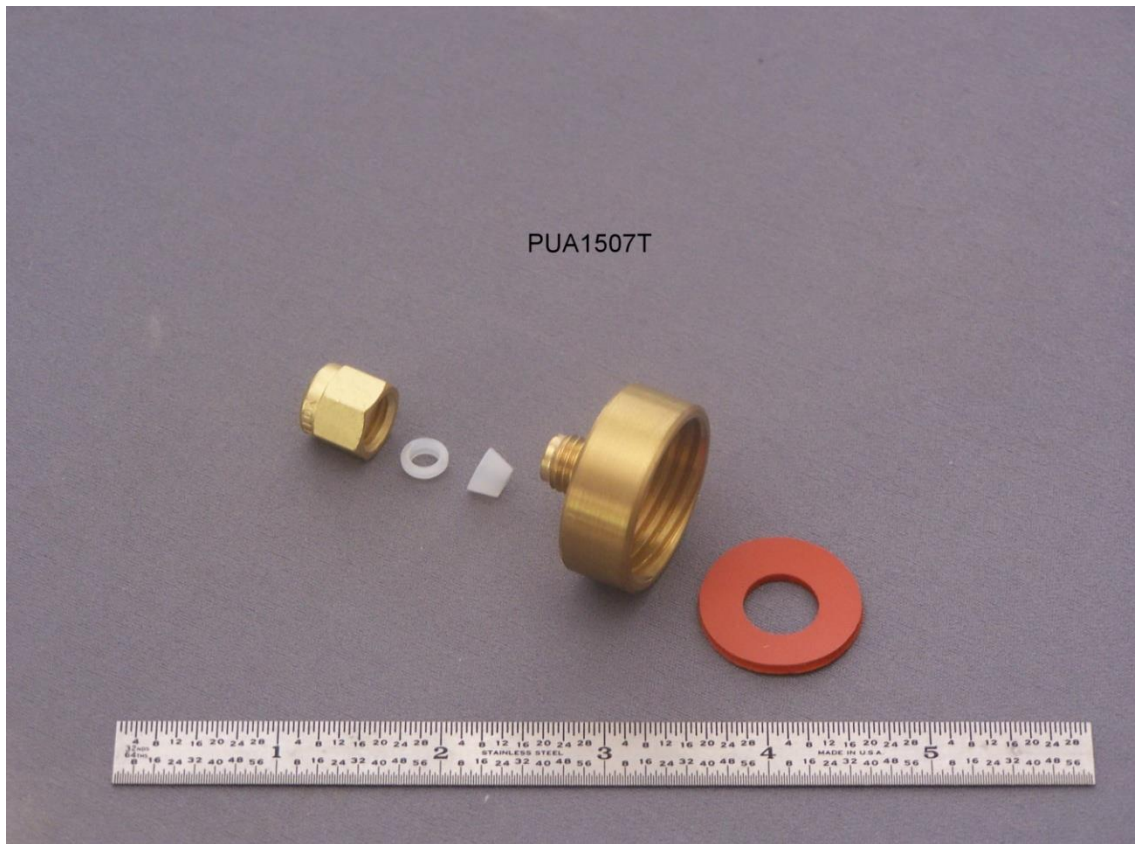

The **Twist Bottle Adapter** provides the means to install a **Sentry** temperature probe into a bottle with a twist-off closure for a test.

To install the probe in the **Twist Bottle Adapter**:

1. Slide the Swagelok nut onto the probe.
2. Slide the flat nylon ferrule onto the probe with the small diameter towards the tip.
3. Slide the graduated nylon ferrule onto the probe with the small end towards the tip.
4. Insert the **Sentry** probe into the **Twist Bottle Adapter** adjusting to the proper height.
5. Tighten the Swagelok nut sufficiently (by hand only) to hold the probe in place.